

# TBS Buffer (20X)

Catalog number: ACC 00580

## LIST OF CONTENTS

Lot number	Item	Concentrate
_____	ACC 00580/0060	60 ml

## STORAGE

TBS buffer (20X) concentrate should be stored refrigerated at 4° C.

Agdia recommends preparing only as much buffer as is needed for one day. Those who store buffers outside of this recommendation are advised to add sodium azide (Sigma S-2002) to 1X liquid buffers at a rate of 0.2 g per liter (0.02%).

## PRECAUTIONS

Prevent direct skin and eye contact with, or ingestion of, product components. Obtain medical attention in case of accidental ingestion of kit components. Always wash hands thoroughly after using this product.

## PREPARING BUFFER

TBS buffer is used in Phytodetek<sup>®</sup> ELISAs for sample preparation, standard dilution and tracer dilution.

Dilute 1 volume of 20X TBS buffer with 19 volumes of distilled water before use.

Buffer concentrate      50 ml

Distilled water          950 ml

To make 1000 ml of TBS buffer, measure the distilled water into a container and add concentrate while stirring.

*Phytodetek<sup>®</sup> is a registered trademark of Agdia, Inc.*

## TECHNICAL SERVICE

If you have any questions about using this product, please contact Agdia, Inc. Monday-Friday by phone (574-264-2014 or 800-622-4342) or by email ([info@agdia.com](mailto:info@agdia.com)).